COLD THAT CARES
evermed

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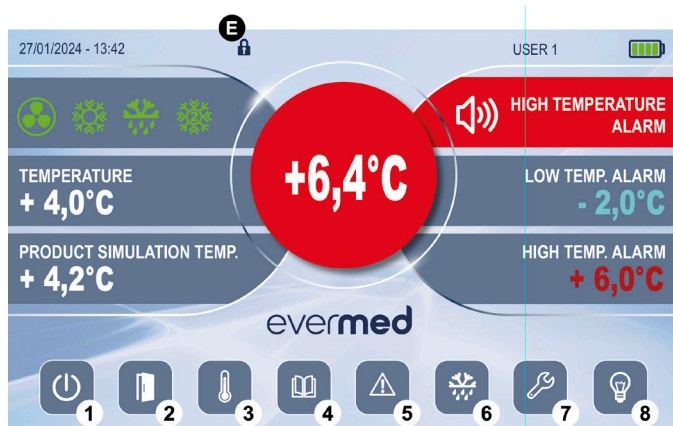
DESCRIPTION OF CONTROLS AND PUSHBUTTONS

DOWNLOAD THE COMPLETE MANUAL

- | | |
|----------------------|---------------------------------------|
| (1) ON/OFF button | (5) Archive/Alarm memory |
| (2) Door button/icon | (6) Start manual defrost (if present) |
| (3) Setpoint button | (7) Menu setting |
| (4) Data menu | (8) Internal lighting (if present) |

START-UP

- When the control panel is powered, immediately after the initial control test, the **ON/OFF** button will appear on a black background.
- press the **ON/OFF** button to switch on the appliance
 - select the language desired by pressing on the corresponding flag icon
 - the alpha-numeric keyboard appears on the display
 - enter the password (**0 by default**) and press ENTER to confirm and turn on the device

STOPPING THE UNIT

- touch the display and enter the password if requested
- press the **ON/OFF** button (1)
- when the pop-up appears, press the confirm button ✓ to turn off the appliance

LOCK/UNLOCK THE DISPLAY

The control panel ADVANCED is equipped with a lock-button protection to prevent tampering by unauthorized personnel that automatically activates within 30 seconds (adjustable) after the last operation carried out on the keyboard itself.

The lock-button condition is indicated by the appearance of the lock symbol **E** at the top of the display. To unlock the control panel, touch the display and enter the password (**0 by default**).

DATE AND TIME SETTING

To set the date and time of the display proceed as follows:

- deactivate any keypad lock
- press the **SETTINGS** button (7) to enter the corresponding menu
- press the **DATE AND TIME** button to open the corresponding pop-up
- select the value to be changed by pressing it (the value is shown blank)
- use arrows to change the value
- validate all changed values by pressing the confirm button ✓

SETTING THE TEMPERATURE

The appliance can be used within the temperature limits specified by the manufacturer.

To set the working temperature to the desired value, proceed as follows:

- from the main menu press the button (3) to open the **SET TEMPERATURE** pop-up
- press the arrow keys to set the desired value of the new temperature
- confirm by pressing the ✓ button or cancel with the X button

The symbol  means that this product should not be treated as a household waste.

To prevent potential negative consequences for the environment and health, be sure this product is correctly disposed of and recycled.

For further information on the disposal and recycling of this product, please contact your Distributor, after Sales Service or the waste Treatment Service.

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